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CHEMICAL SPILLS RESPONSE KIT USAGE INFORMATION

1- Personal protective equipment is worn.

2- Warning sign is put.

3- The chemical spill is surrounded by an absorbent sausage, the spread of the waste is prevented.

4- An absorbent pad is placed on the chemical waste, chemical waste is expected to be fully absorbed.

5- Remaining residues in the area, the dust absorbent is poured and it is solidified and collected with the help of a dustpan and a scraper.

6- Solidified residue wastes, used pads, sausages, are put into a hazardous waste bag.

7- Spillage area is cleaned with a cleaning sponge.

8- Cleaning is terminated by cleaning the area thoroughly with cleaning clothes.

9- All equipment including personal protective equipment used in cleaning the materials used are put into the hazardous waste bag and fastened with a plastic clamp, sent to the hazardous waste storage.

- In case of emergency call **1122** immediately.
- If anything you need, contact the lab administrator.
- If you have questions about the spill kit, don't hesitate to mail isg@ku.edu.tr.